

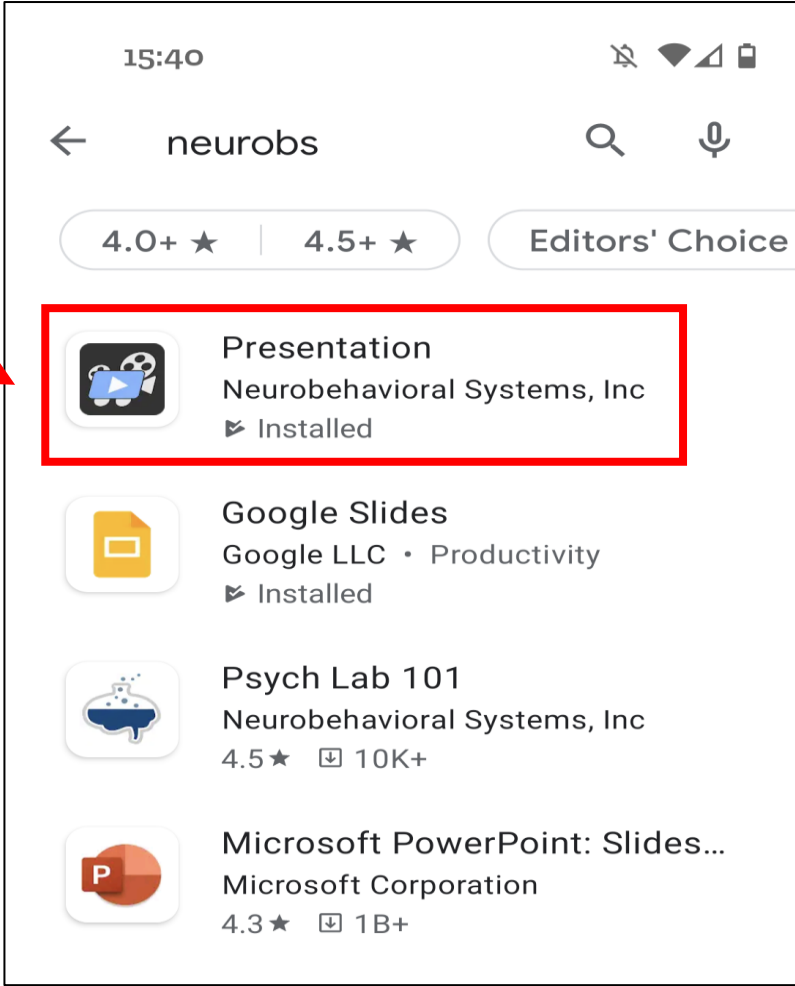


Step 1: Go to google play store/ app store, search for **Neurobehavioral Systems** and select **Presentation**

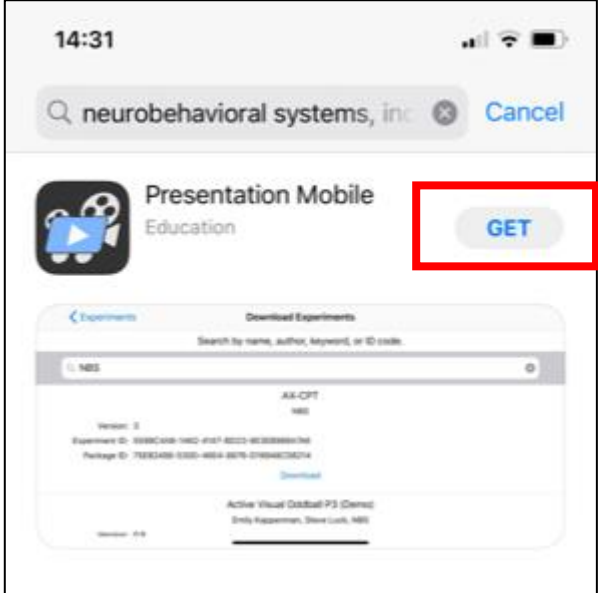


Step 2: Press **Install / Get**

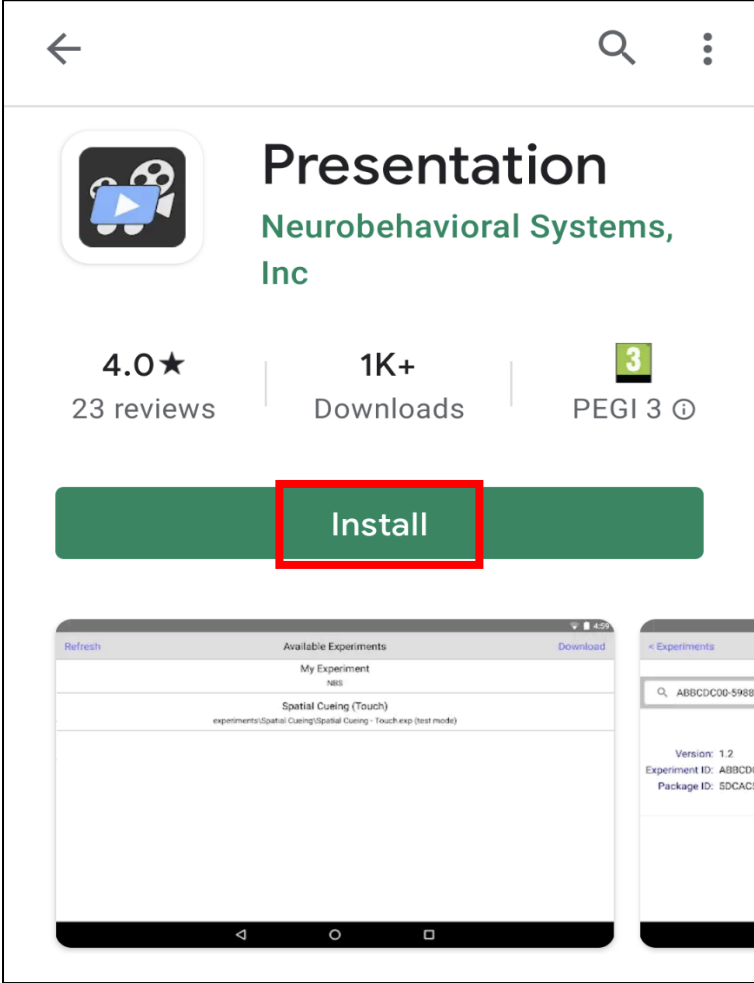


Step 3: Open the application. Keep in mind under *Installed Experiments* there is nothing yet because we haven't downloaded anything.
Now select **Download**




Step 4: Type iPRESSmobile on the search bar and press Search

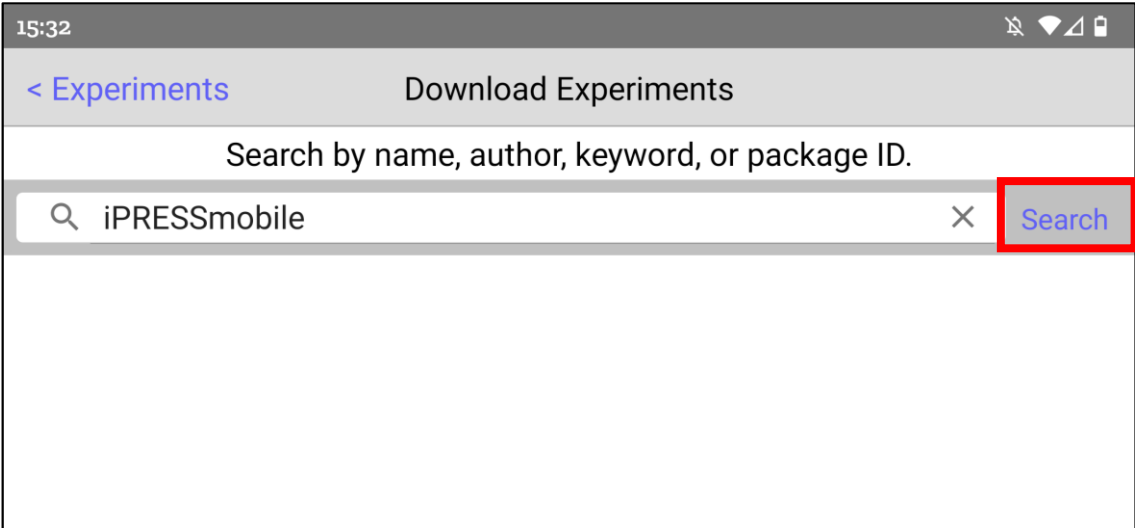


The search results will show **WMU Training Practice (mobile)**

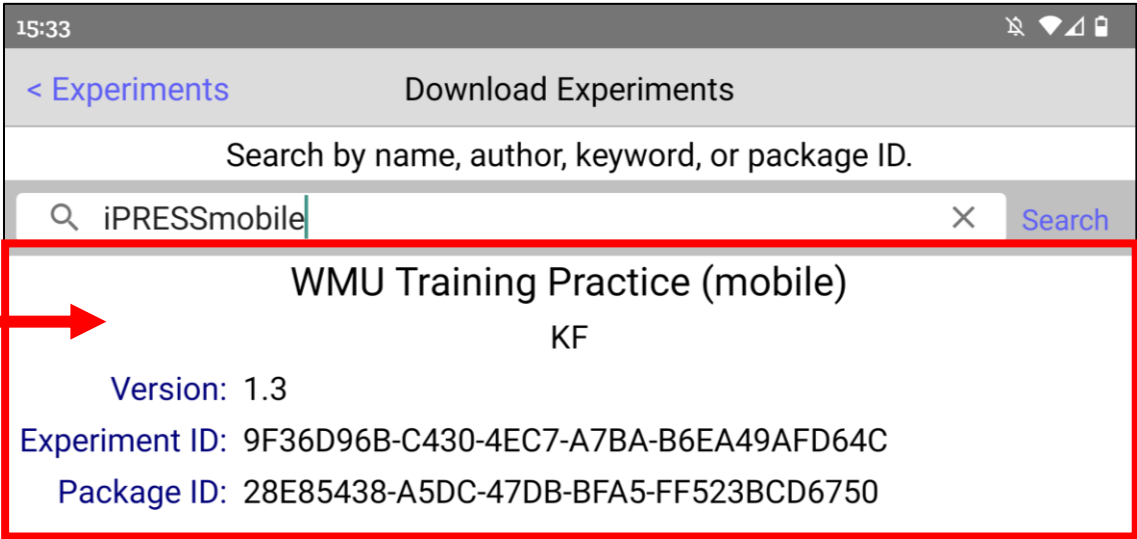

Step 5: Scroll down and press **Download** at the bottom

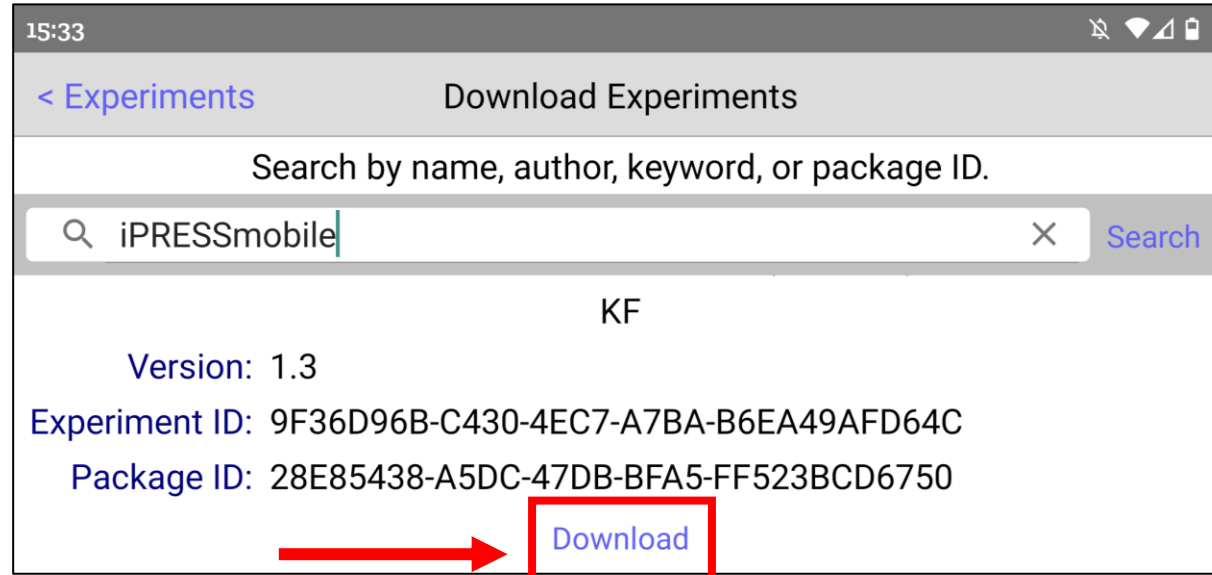

Now the WMU Training Practice experiment will appear under ***Installed Experiments***

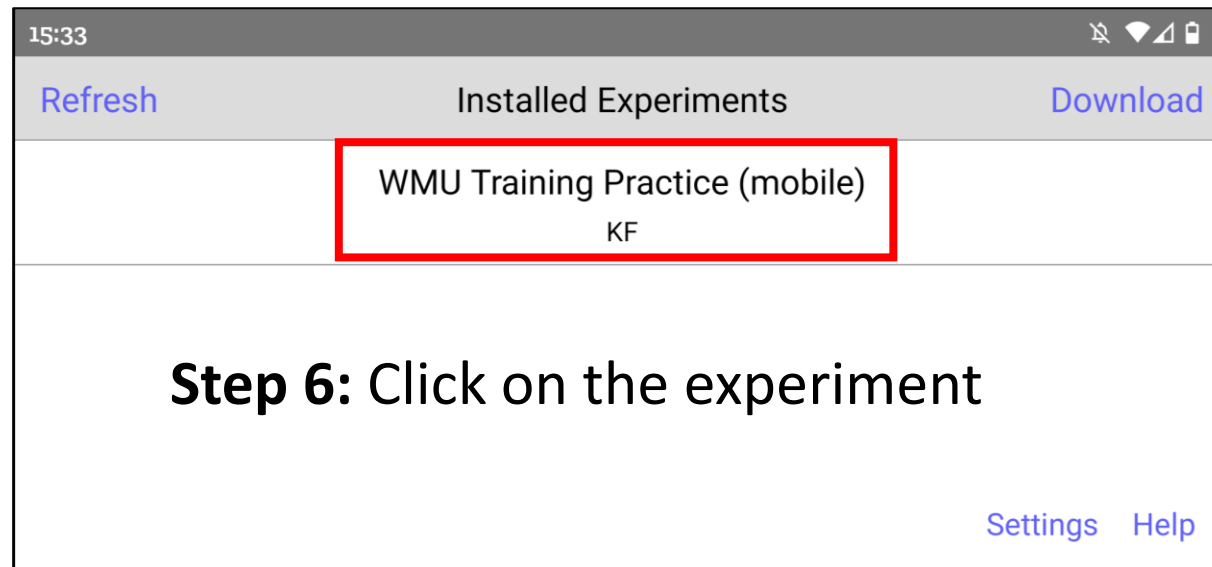

Step 7: Write your unique **Subject ID** (I will let you know what that is)

15:33

< Experiments WMU Training Practice (mobile) Info

Scenario:
MatrixPractice1.sce
MatrixPractice2.sce
Matrix_withFeedback.sce

Subject ID: test

Start

(During experiment, touch upper-left and lower-right corners simultaneously to quit.)

In the example I inserted **“test”** as the ID

Step 8: Press **Start** to begin the experiment

15:34

< Experiments WMU Training Practice (mobile) Info

Scenario:
MatrixPractice1.sce
MatrixPractice2.sce
Matrix_withFeedback.sce

Subject ID: test

Start

(During experiment, touch upper-left and lower-right corners simultaneously to quit.)